Technical Sheet

N° 401476 BI -0 02/2010 pag. 1 / 2

MODIFIED LAURYL SULFATE TRYPTOSE (mLST) BROTH BASE

ENTEROBACTER SAKAZAKII ISOLATION AGAR (ESIA)

Enrichment selective broth and selective chormogenic plating medium for the detection of Cronobacter (Enterobacter) sakazakii in milk and milk products.

TYPICAL FORMULATIONS

mLST Broth Base –			
POWDERED MEDIUM (g/L)		ESIA - POWDERED EMDIUM AND READY TO USE PLATES(G/L)	
PEPTONE*	20.00	PEPTONE	7.00
LACTOSE	5.00	YEAST EXTRACT	3.00
POTASSIUM PHOSPHATE MONOBASIC	2.75	SODIUM CHLORIDE	5.00
POTASSIUM PHOSPHATE BIBASIC	2.75	SODIUM DESOXYCHOLATE	0.60
SODIUM LAURIL SULFATE	0.10	CRYSTAL VIOLET	0.002
SODIUM CHLORIDE	34.00	X-α-GLUCOSIDE	0.150
* Enzymatic digest of animal and plant tissue		A GAR	15.00

VANCOMYCIN ANTIMICROBIC SUPPLEMENT (x vial)

VANCOMYCIN 25 mg

mI CT	DDOTH	DE 4 DV TO		(~/L)
misi	BROTH -	READY TO	USE TUE	RES (0/L)

Peptone*	20.00
LACTOSE	5.00
POTASSIUM PHOSPHATE MONOBASIC	2.75
POTASSIUM PHOSPHATE BIBASIC	2.75
SODIUM LAURIL SULFATE	0.10
SODIUM CHLORIDE	34.00
VANCOMYCIN	0.01
* Enzymatic discost of animal and plant tipous	

Enzymatic digest of animal and plant tissue

METHOD OF PREPARATION

Modified Lauryl Sulfate Tryptose (mLST) Broth Base

Suspend 64.6 g in 1000 ml of cold distilled water. Heat to dissolve and sterilize by autoclaving at 121° C for 15 minutes. Cool to room temperature and add 2 ml of the vial contents of Vancomycin Antimicrobic Supplement (REF 4240057), reconstituted with 5 ml of sterile distilled water (the reconstituted supplement can be stored at -20 °C for 1 month). Mix well and distribute into sterile tubes (10 ml/tube) final pH 6.8 \pm 0.2

Enterobacter Sakazakii Isolation Agar (ESIA)

Suspend 30.8 g in 1000 ml of cold distilled water. Heat to boiling stirring constantly; sterilize by autoclaving at 121 °C for 15 minutes. Cool to 45-50 °C and distribute into sterile petri dishes. final pH 7.0 ± 0.2

TECHNIQUE

- To prepare the primary dilution, add x g of the test sample to 9 times x ml of Buffered Peptone Water (REF 401278).
- Allow dry samples to disperse in the liquid without stirring. If a sample has not been dissolved completely after 30 min, than mix it gently with the medium.
- Incubate the inoculated Buffered Peptone Water at 37 °C ± 1 °C for 18 h ± 2 h.
- After incubation of the inoculated pre-enrichment medium, transfer 0,1 ml of the obtained culture into 10 ml of mLST/vancomycin medium.
- Incubate at 44 $^{\circ}$ C ± 0,5 $^{\circ}$ C for 24 h ± 2 h.

Biolife Italiana Srl

Technical Sheet

N° 401476 BI -0 02/2010 pag. 2 / 2

- It is recommended to use either a water bath (6.3) or a forced-air incubator to ensure that the maximum temperature (44,5 °C) is not exceeded.
- After incubation of the inoculated mLST Broth with vancomycin streak a loopful (ca. 10 μl) onto the surface of the ESIA plate.
- Incubate the plate at 44 °C ± 1 °C for 24 h ± 2 h.
- After incubation, examine the chromogenic plate for the presence of typical colonies of presumptive Cronobacter (Enterobacter) sakazakii.
- Typical colonies are small to medium sized (1 mm to 3 mm) green to blue-green colonies. Non-typical
 colonies are often slightly transparent and violet coloured.
- Confirm the colonies with the biochemical tests recommended by ISO/TS 22964.

User Quality Control (mLST Broth)

Productivity: *C.sakazakii* good growth Selectivity: *S.aureus* ATCC 25923: no growth Incubation at 44 $^{\circ}$ C \pm 0,5 $^{\circ}$ C for 24 h \pm 2 h

User Quality Control (ESIA)

Productivity: C.sakazakii good growth, blue colonies

Specificity: Enterobacter gergovieae: good growth, purple colonies

Selectivity: S.aureus ATCC 25923: no growth Incubation at 44 $^{\circ}$ C $^{\pm}$ 0,5 $^{\circ}$ C for 24 h $^{\pm}$ 2 h

STORAGE

mLST Broth (powdered medium): 10-30 °C

ESIA (powdered medium): 2-8 °C Ready to use tubes and plates: 2-8 °C

LIMITS

Some coliform bacteria grow on ESIA plates with violet colonies.

REFERENCES

•ISO/TS 22694:2006 Milk and milk products - Detection of Enterobacter sakazakii

PΑ	CK	٩GI	NG

4014762	Modified Lauryl Sulfate Tryptose (mLST) Broth Base	500 g (7,8 L)
4014782	Enterobacter Sakazakii Isolation Agar (ESIA)	500 g (16,2 L)

SUPPLEMENT

4240057	Vancomycin Antimicrobic Supplement	10 VIALS
---------	------------------------------------	----------

5514762	Modified Lauryl Sul fate Tryptose (mLST) Broth	20x10 ml (ready to use tubes)
541478	Enterobacter Sakazakii Isolation Agar (ESIA)	20 ready to use plates