

## VANCOMYCIN ANTIMICROBIC SUPPLEMENT

**VIAL CONTENTS**

Vancomycin.....25 mg

**DIRECTIONS**

Prepare 1000 ml Modified Lauryl Sulfate Tryptose (mLST) Broth Base (REF 4014762) autoclaved at 121 °C and cooled to room temperature. Aseptically reconstitute the contents of one vial of Vancomycin Antimicrobial Supplement with 5 ml of sterile distilled water. Add 2 ml of reconstituted supplement to 1000 ml of mLST Broth. Mix well and aseptically distribute into sterile tubes (10 ml/tube).

The remaining unused supplement can be stored at -20 °C for 1 month.

**INTENDED USE**

For the preparation of the enrichment selective broth for the detection of *Cronobacter (Enterobacter) sakazakii* in milk and milk products.

For the details of the procedure see the technical sheet of Modified Lauryl Sulfate Tryptose (mLST) Broth Base (REF 4014762).

**STORAGE**

Store at 2-8 ° - When stored as directed the supplement remains stable until the expiry date shown on the label. Do not use beyond stated expiry date.

**PRECAUTIONS AND WARNING**

For Laboratory use only. The supplement should be used only by adequately trained personnel with knowledge of microbiological techniques in the laboratory.

Consult the material safety data sheet before the use.

**PACKAGING**

4240057 Vancomycin Antimicrobial Supplement

10 vials